



Precision weighing instruments since 1880

1401 17th Street, Suite 750
Denver, CO 80202
800.321.1135
303.431.7255
303.423.4831 FAX
www.denverinstrumentusa.com

Application Note

Setting Up Dot Matrix Printer with 200 Series Meter

Printer settings for Denver dot matrix printer 902579.1 and Denver 200 Series meter with default serial settings.

To change the DIP Switch:

1. Remove the clear cover.
2. Look inside the printer in the direction of the keys.
3. The DIP switch is blue with 6 white switches.
4. Cycle power after changing any DIP switches.

NOTE: "ON" means away from the paper. "OFF" means towards the paper and nearest to the number on the DIP switch.

NEW 200 Series meters (serial numbers beginning with letter E):

Switch 1 – ON
Switch 2 – OFF
Switch 3 – OFF
Switch 4 – ON
Switch 5 – ON
Switch 6 – OFF