DLT Series

Operating Instructions | Betriebsanleitung | Mode d'emploi | Manual de instrucciones | Istruzioni per l'uso | 取扱説明書 | 操作手册

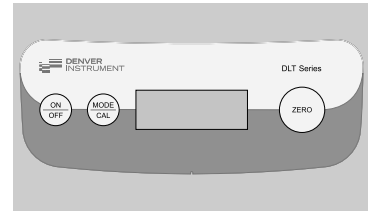
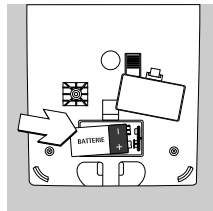
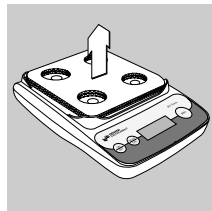
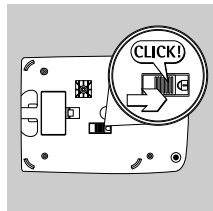
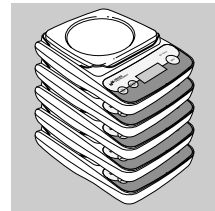
Electronic Precision Balances | Elektronische Präzisionswaagen | Balances électroniques d'analyse et de précision | Balanzas electrónicas de precisión | Bilance elettronica di precisione | 電子上皿天びんイーコンシリーズ | 精密電子天平

Europa, Asia und Australia:
Robert-Bosch-Breite 10
37079 Goettingen, Germany

North and South America:
1401 17th St. Suite 750
Denver, CO 80202, U.S.

www.denverinstrument.com

GB/U.S.

Warnings and Safety Precautions

- To prevent damage to the equipment, please read these operating instructions carefully before using your balance.
- Do not use this equipment in hazardous areas.
- Exposure to excessive electromagnetic interference can cause the readout value to change. Once the disturbance has ceased, the instrument can be used again in accordance with its intended purpose.
- Do not open the balance housing as this may void the manufacturer's warranty if the seal is damaged.

Storing the Balance

- Do not stack more than 5 balances on top of one another for storage.

Transportation Locking Device 1

- Unlock the balance as follows: Move the switch in the direction of the arrow

Transportation Locking Device 2

- Remove the cardboard between the balance and the weighing pan
- Reposition the weighing pan

AC Power Connection

- Optional AC adapter: Please contact your local dealer

Installing the Battery

- Use only commercially available 9-volt batteries (included in the standard equipment supplied) or a rechargeable battery. Inserting the wrong type of battery can result in an explosion.
- If you use a rechargeable battery, recharge it only with an external battery charger.
- Open the battery compartment: remove the compartment cover
- Install the 9-volt battery or rechargeable battery in the battery compartment
- Make sure the polarity is correct.
- Close the battery compartment: press down on the compartment cover to snap it into place

Operation: Basic Weighing Function | Functions of the Keys

To switch on or off: → (ON/OFF) → → (ON/OFF) → Press and hold for more than 2 seconds

To zero: → (ZERO) → → (ZERO) →

To toggle weight unit: → (MODE) → → (MODE) → → (MODE) → → etc.: (MODE) →

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GB/U.S.

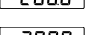
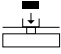
Automatic Shut-off

- Battery operation: automatic shut-off after 6 minutes
Example:
If the weight readout remains unchanged and no keys are pressed for at least 4 minutes, the display starts flashing. After another 2 minutes, the balance will shut off automatically unless a key is pressed.

Calibration/Span Adjustment

- Purpose: Calibrates and adjusts the span of the balance after it is initially powered on for optimal accuracy
- Optional calibration weight: please contact your local dealer

Calibration/Span Adjustment

- Zero the balance:**  (ZERO) → 
- Start calibration by pressing and holding (CAL) for 5 seconds** → 
- Place the required calibration weight on the balance** (not included in std. equipment supplied): 
- Complete calibration:** Remove the calibration weight when "g" is displayed next to the weight. 

Error Codes

Display/Problem	Cause	Solution
No segments appear on the display	Battery is drained	Replace battery, recharge battery using external charger
-H-	The load exceeds the balance capacity	Unload the weighing pan
-L-, Err56	Weighing pan not in place Preload during power-up too high	Place the weighing pan correctly on the balance Unload the pan and switch on the balance again The weighing pan may not touch any surrounding items
Err2	Calibration condition not met, e.g.: - Balance not zeroed! - The weighing pan is loaded	Unload balance and press (ZERO) Calibrate only when zero is displayed
Max. weighing capacity is less than indicated under "Specifications"	The balance was switched on without the weighing pan in place	Place the weighing pan on the balance and press ON/OFF
The weight readout is obviously wrong	The balance was not calibrated/adjusted Balance not zeroed	Calibrate/adjust the balance Zero the balance

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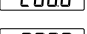
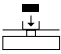
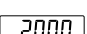
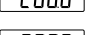
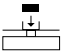
Specifications

Model	DLT-411	DLT-211	DLT-4100	DLT-2100
Weighing capacity/readability	g 410/0.1	210/0.1	4,100/1	2,100/1
Linearity	≤±g 0.2	0.2	2	2
Operating temperature range	+10... +30°C (273 ... 303 K, 50° ... 86 °F)			
Calibration weight	g 200	200	2000	2000
Net weight Pan size	440 g 135×140 mm			
Power source Voltage Frequency	Optional power supply 230 V/115 V, +15%...-20%			
Input voltage range	7 V to 20 V; power consumption (average)			
Hours of operation w/9-volt battery (alkaline) - Rechargeable battery fully charged (NiMH)	approx. 30 h 8 h (avg.)			

CE, CSA, UL approval depending on area where applicable

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